



## ChIP Validated Sin3a Antibody with Positive Primer Set

Catalogue no: 900020

Chromatrap®'s ChIP Validated Sin3a Antibody and Positive Primer Set provides a complete set of tools to assist with a successful ChIP assay. Including: Sin3a antibody, control sheep IgG, and a positive primer set. The ChIP Validated Sin3a Antibody and Positive Primer Set is not suitable for use with non-human species.

### Background:

Sin3a (Switch independent 3 homologue) is a transcriptional repressor and protein-encoding gene. It interacts with co-repressors and DNA-binding proteins such as REST, MXI1 and MXD1-MAX that allow Sin3a to bind to DNA and repress transcriptional activity. It acts interdependently with OGT (O-linked N-acetylglucosamine) to repress transcription alongside with histone deacetylation.

A sheep IgG is included in this Antibody Primer Set as a negative control for the ChIP experiment.

The Sin3a positive primer set recognises the hypermethylated promoter of a gene that binds to Sin3a, resulting in the repression of transcription.


### Suggested Usage:

Component	Suggested Dilution	Figure
Sin3a	2:1 (antibody: chromatin)	1
Sheep IgG	2:1 (antibody: chromatin)	1
Positive Primer Set	Dilute from 4µM (provided) to 1µM working concentration	

Please note: Optimal dilutions should be determined by the user. These volumes are stated as guidelines only.

### Fig 1. Sin3a ChIP qPCR

Chromatin immunoprecipitation (ChIP) assays were performed using the Chromatrap® standard ChIP spin column sonication kit for qPCR (Cat no. 500071) with 1ug of chromatin from MCF7 cells and 2ug of Anti- Sin3a antibody. qPCR was used to analyse the enrichment of Sin3a onto the positive gene locus compared to the negative gene locus.

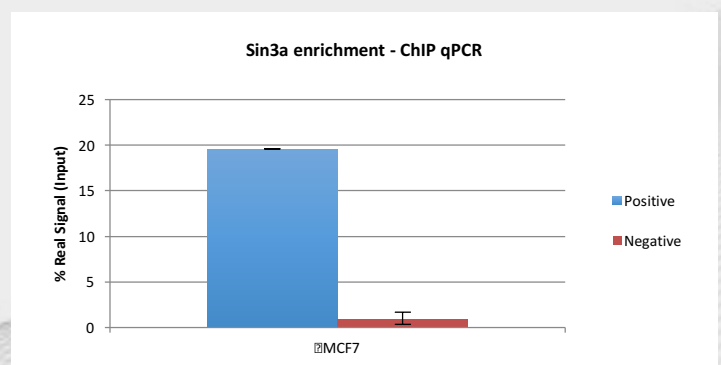

## Advancements in Epigenetics

\*This product is for research use only. There is a possibility that results may vary between antibody lots.



[www.chromatrap.com](http://www.chromatrap.com)

Clywedog Rd South  
Wrexham Industrial Estate  
Wrexham LL13 9XS, United Kingdom

Tel: +44 (0) 1978 666239/40  
Email: [support@chromatrap.com](mailto:support@chromatrap.com)

**Applications:** CHIP

**Concentration:** 1mg/ml

**Size:** 50 $\mu$ l

**Specificity:** Human

**Source:** Sheep

**Type:** Monoclonal

**Purification:** Protein G (affinity purified)

**Storage Conditions:** The Sin3a antibody and sheep IgG should be stored at +4°C. The positive primer set should be stored at -20°C (*Avoid multiple freeze/thaw cycles as this may denature the antibody and degrade the primer sets*)

**Advancements in Epigenetics**

\*This product is for research use only. There is a possibility that results may vary between antibody lots.

Company registered in England no 2047818 Registered office: 7 Regis Place, Bergen Way, Kings Lynn, PE30 2JN United Kingdom